


HIGH RESOLUTION PEPTIDE MAPPING


Tryptic Maps of Transferrin

HAISIL 100

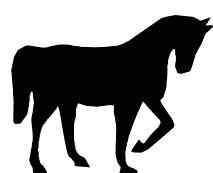

Wide Format Variety:

- ✓ 2cm - 25cm Length
- ✓ 170 μ m - 20mm I.D.
- ✓ Analytical Guard Cartridges
- ✓ Prep Guard Cartridges

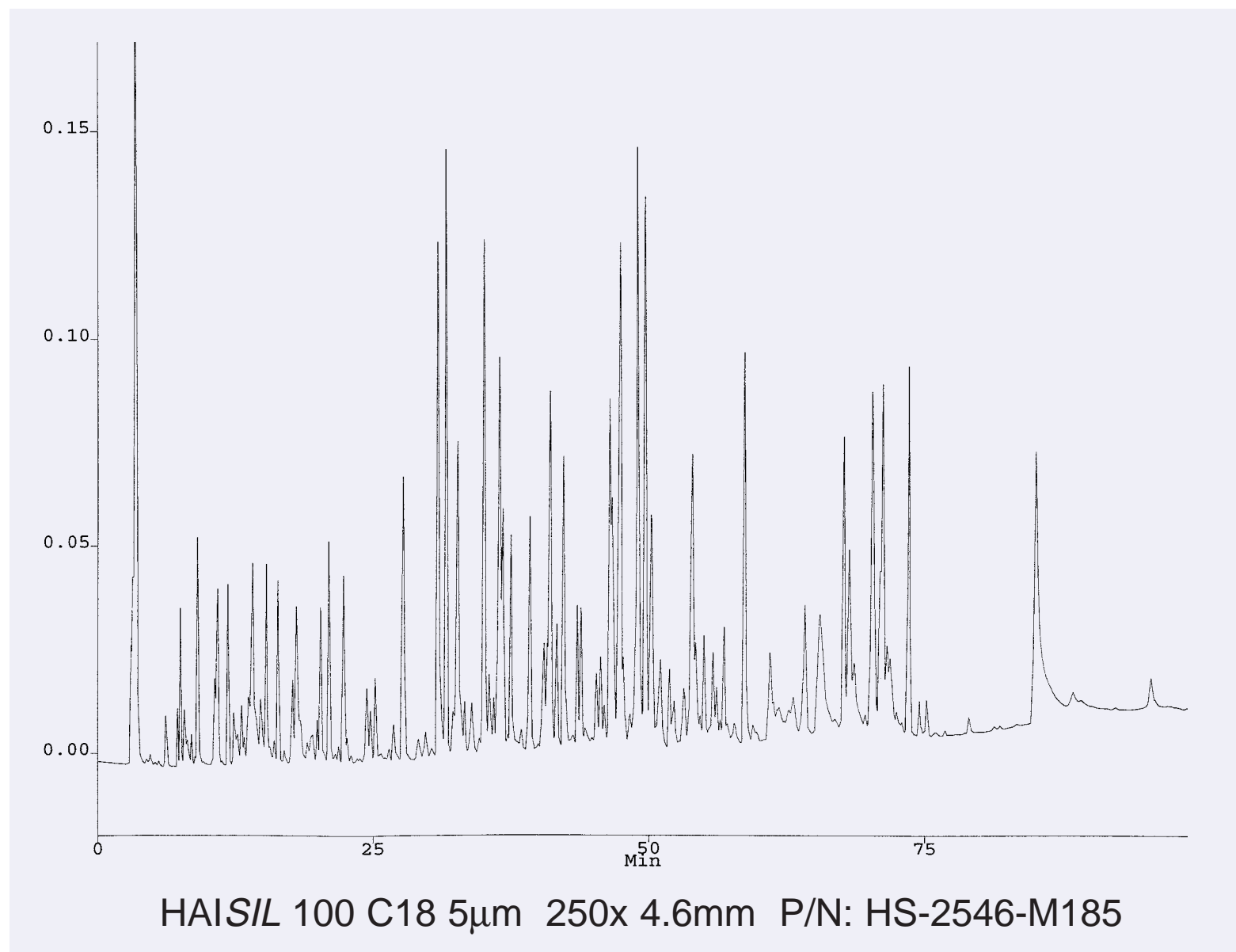

HAISIL 300



Chromatographic Conditions:

Mobile Phase A:
50mM phosphoric acid pH 2.8

Mobile Phase B:
60% MeCN containing
50mM phosphoric acid pH 2.8

Gradient:
0%B for 5min
0-40%B in 120min
40-60%B in 10min

Flow Rates:
900 μ L/min on 4.6mm I.D. column
200 μ L/min on 2.1mm I.D. column

